SM series microscope

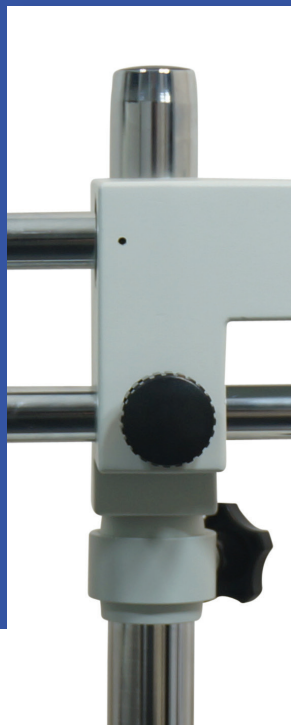
ASSEMBLY MANUAL

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MONTAGE HANDLEIDING

THANKS FOR YOUR PURCHASE

Thank you for purchasing this microscope. We hope you enjoy using it! This manual contains assembly instructions for the SM-4TP and SM-4T microscopes. The assembly process is similar for the models with a single arm.

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01 SET UP THE BASE

The first step of the assembly is setting up the base. There is a small screw present in the base which should not protrude the vertical rod. Make sure that the small locking screw is fully unscrewed before inserting the vertical rod.

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02 TIGHTEN THE VERTICAL PART OF THE STAND

The next step is tightening the vertical part of the stand (1). Lock the thread of the vertical rod into the base. Make sure you tighten the rod firmly to ensure a stable stand. After you tightened the rod firmly, tighten the locking screw again (2). The ring, see (3), is meant to hold the horizontal arm up while you adjust it.

Make sure the ring is securely fastened before you execute the next step.

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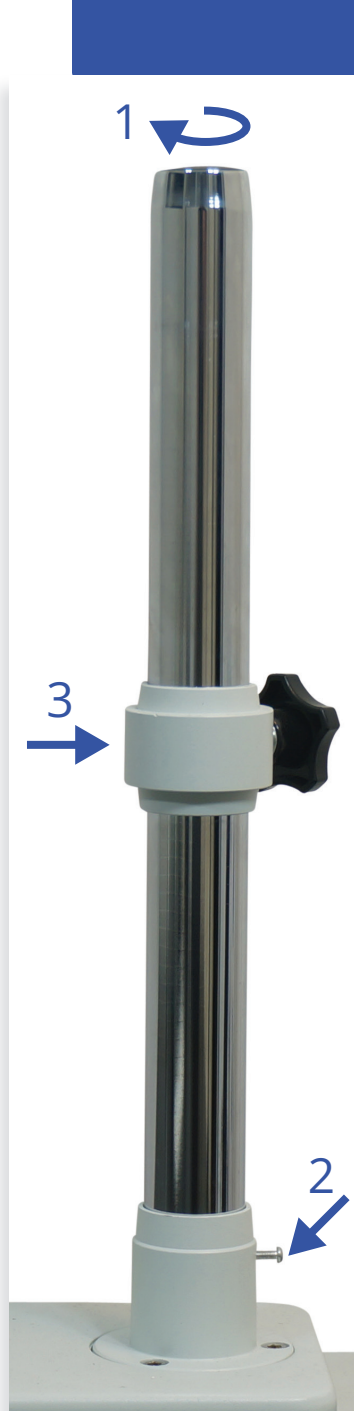
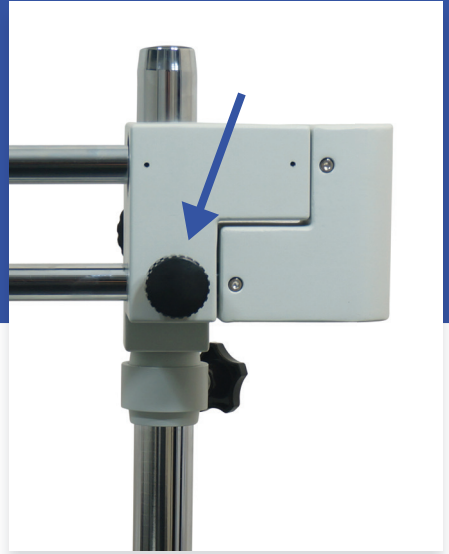
03 ATTACH THE ARM

Lower the arm of the microscope onto the vertical rod. The arm has two screw holes to secure the arm vertically and horizontally during use.

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04 ATTACH THE LAST PART OF THE STAND

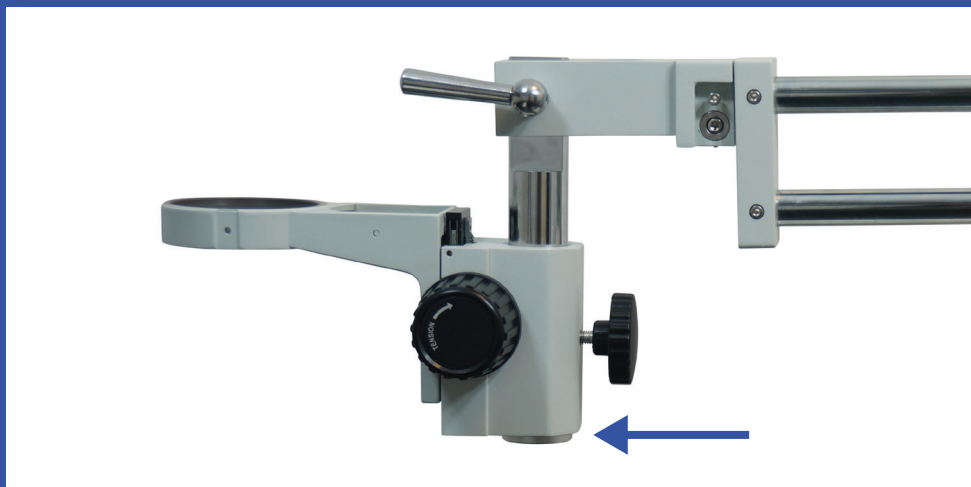
The next step is attaching the last part of the stand. The arrow in the image indicates the place of the two hex screws. Unscrew these first to push the last part of the microscope over the end of the arm. The small screw prevents the part from falling off the stand, the large screw prevents rotation.

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05 MOUNT THE MICROSCOPE HEAD HOLDER IN PLACE

The arrow in the image above indicates the place of the screw cap. It needs to be removed to mount the microscope head holder in place. Unscrew the screw cap anticlockwise and remove it. After that, slide the head over the rod and screw the cap back on. This prevents the head from falling when the setscrew is loosened.

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06 ATTACH THE MICROSCOPE HEAD

The microscope head can now be mounted in the holder. The thumbscrew, indicated by the arrow in the image, can be tightened in 3 places to fixate the head. Use one of the three holes to tighten the head.

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07 THE FINAL ASSEMBLY STEP

Finally, the rubber eyepieces can be placed on the oculars. Small black screws are present in the oculars. These need to be loosened first and then fixated again when the eyepieces are placed. The oculars can be focused individually so that the people with glasses can adjust each eye separately. Before use, remove the protective cap on the bottom of the microscope head.

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08

HOW TO FOCUS

First adjust the height of the microscope head to focus the lenses roughly, then use the button (at the head) to finetune the focus of the lenses. For more freedom of movement, use the large screw on the back of the microscope head holder.

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ACCESSORIES (SOLD SEPARATELY)

Microscope camera

A camera can be added to the SM series microscopes. Please note that you can only use a camera in combination with a mount adapter, which is also sold separately.

(C-) mount adapter

If you would like to use a camera on your microscope, you need an adapter to attach the camera to the microscope.

Barlow lens

Use a Barlow lens to increase the distance between the microscope head and the object you want to inspect. This way, you can increase the height of your workspace underneath the microscope.

LED ring

The microscope head can cast a shadow on the object that you are inspecting. In addition, lighting might not always be ideal. To get great lighting on your object, attach a LED ring to the head of the microscope.

Premium rubber eyepieces

For more comfort while looking through the oculars, use premium rubber eyepieces. The eyepieces also reduce light incidence.

Oculars

Sets of oculars with different magnification levels can be purchased separately. For example: oculars with a magnification of 15x or 20x instead of the 10x magnifying oculars that are standard included with SM series microscopes.

For more information or the latest updates,
please check elishop.eu/microscope

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